



※ Prior to operate the Maelstrom 4810, read carefully the User Manual is a must.

STEP 1 Power on

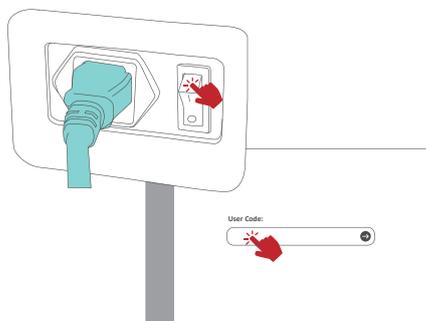


Power on the system and login by default User Code: 333.

STEP 2 Tap Tip icon

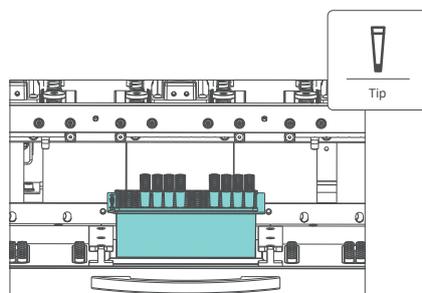


Place one tip box, then tap the "Tip" icon.
PS: Only one tip box at a time

STEP 3 Mount Tips

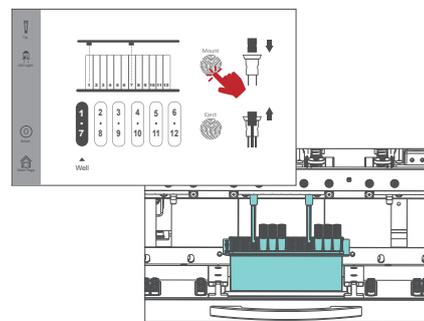


Choose the tip position then tap "Mount".

STEP 4 Place Samples

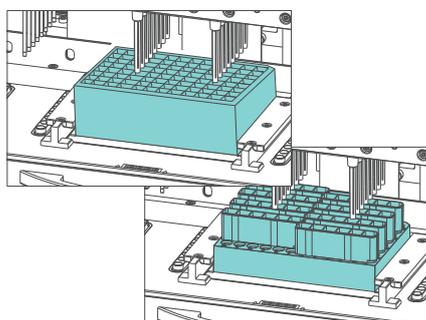


Replace with pretreated Auto Plate/Tube.

STEP 5 Start the Run

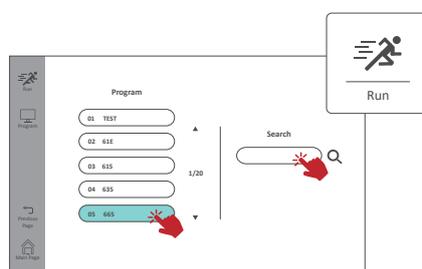


Press on "Run" and double tap the desired program.

STEP 6 Run to Finish

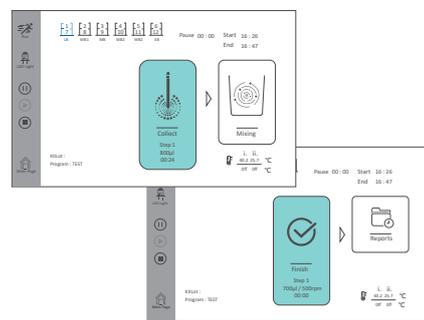


Follow the on-screen instruction to start the extraction.

STEP 7 Eject Tips

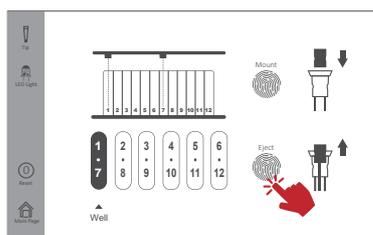


Tap "Tip" → "Eject" to eject the tips if choosing Auto Tube. Withdraw the finished plate.

STEP 8 Decontamination

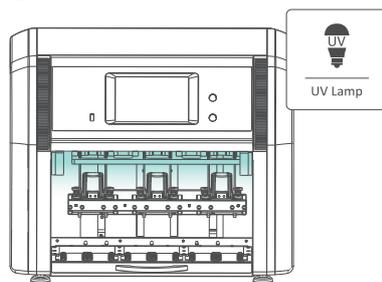

Turn on "UV Lamp" for 30 mins after extraction.

Maelstrom 4810

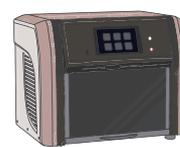

Taiwan Advanced Nanotech Inc. www.tanbead.com