




# Switch 8

## Import Programs Instruction

[www.tanbead.com](http://www.tanbead.com)

Step 1. Login with 222 as Supervisor to access Edit in the main menu

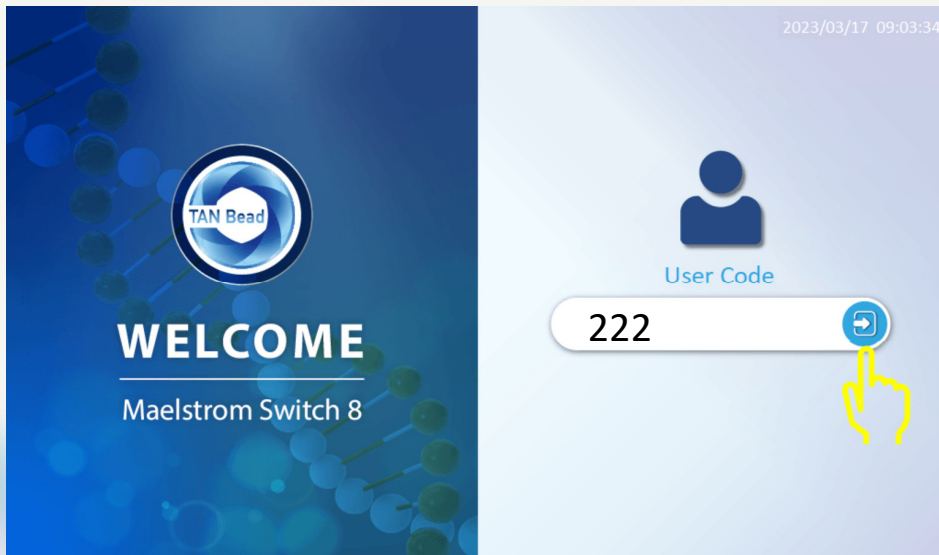

TANBEAD

Step 2.


Insert the USB flash drive with the program files, screen will display the following message



TANBEAD

Step 3.


Choose the Program number for the new program position, then press Import

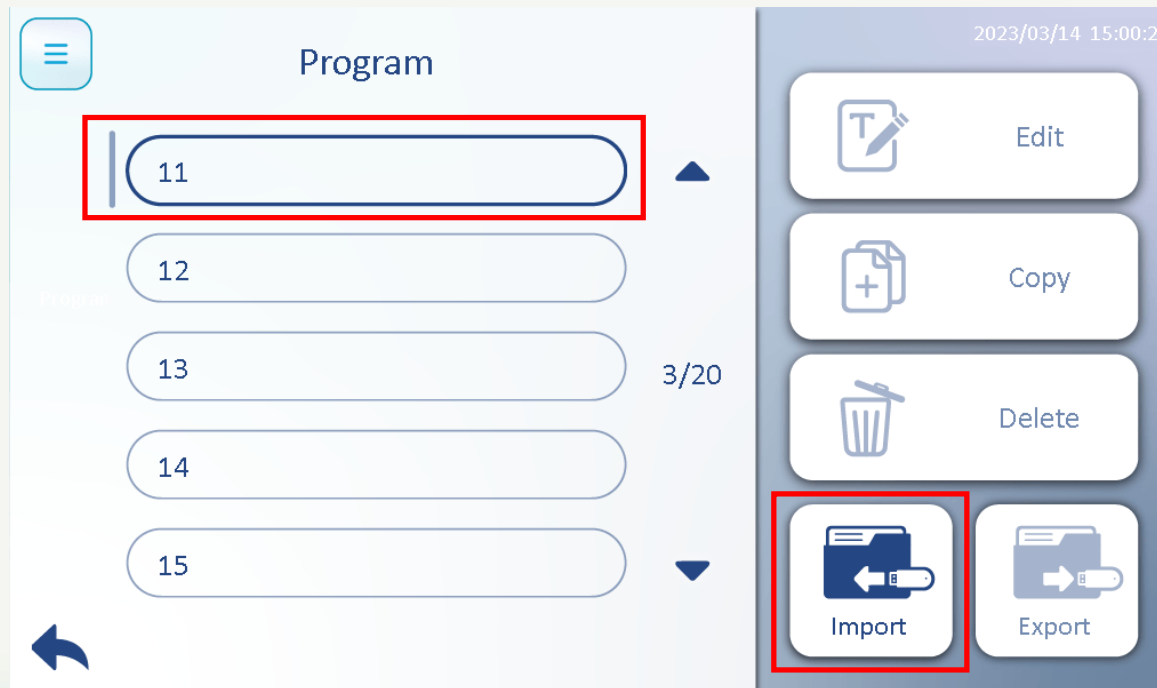


TANBEAD

Step 4.

Select the program file from the USB flash drive, then press “Open” to load the new program.

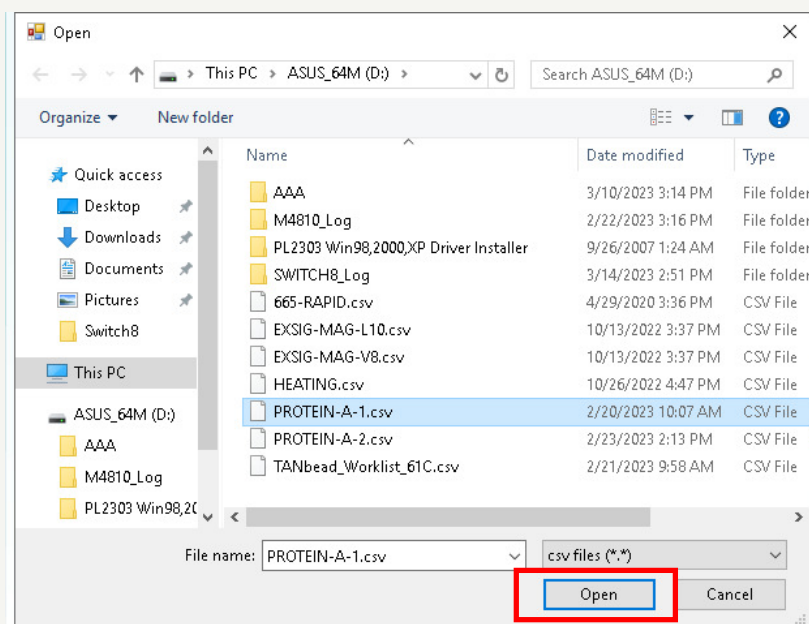


05 613

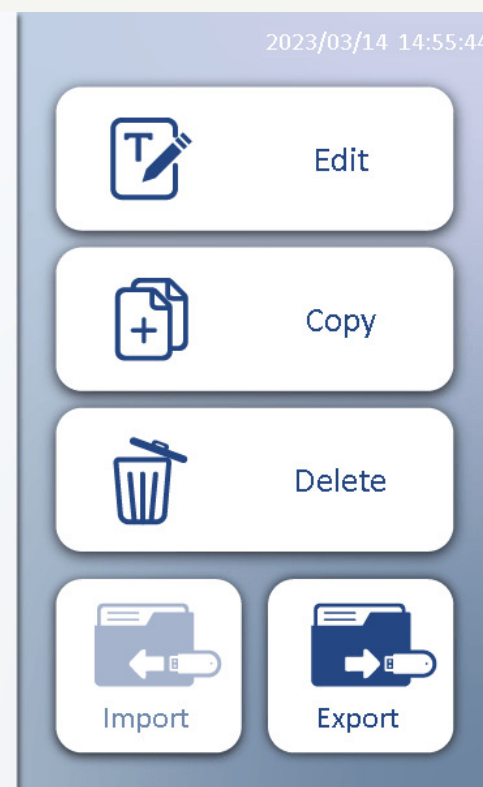

TANBEAD

Step 5.

A Success message will pop up, then you can find the new Program has been imported to the designate program position

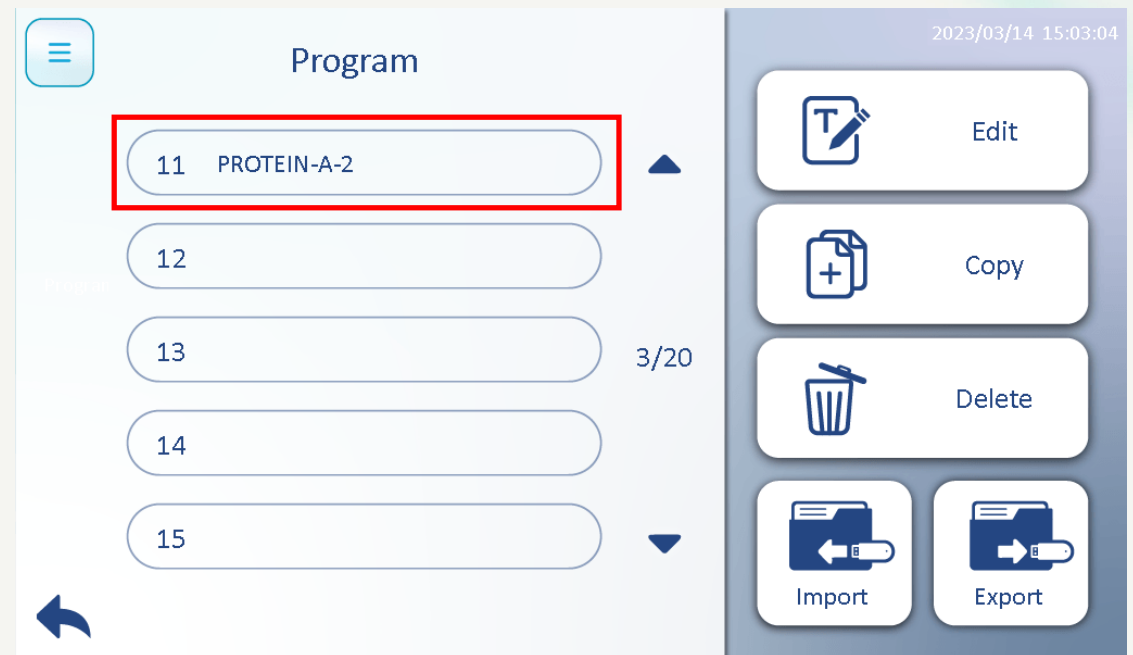

TANBEAD