



TANBead® Nucleic Acid Extraction Kit

Viral Auto Tube

(For use with the SLA-16 / 32 / E13200 series)

RUO

635S46

(For Research Use Only) V1

1. Intended Use

This product is designed for isolating nucleic acid from various samples, which can be performed by using TANBead® Nucleic Acid Extractor and is intended for research use only.

2. Purpose

TANBead® Nucleic Acid Extraction Kit (635S46) is suitable for isolating a variety of viral nucleic acids such as Enterovirus, Japanese encephalitis virus, Dengue virus, Avian Influenza virus, and Epstein-Barr virus. Serum specimens are processed through a series of automatic extraction steps and finally the high-quality nucleic acids can be applied directly to the following qualitative and quantitative assays. With high sensitivity, this reagent kit can be applied for clinical diagnostics and research.

3. The basic principle

The silicon dioxide layer coated on the magnetic beads can adsorb the negatively charged molecules to purify nucleic acids from samples.

4. Specification

Starting Materials	200 µL serum or PBS suspension
Elution Volume	70~100 µL

5. Component Supplied with the Kit

Auto Tube	8 trays	Auto Tube with reagent buffers
Proteinase K	1.0 mL	Proteinase K
Elution Buffer	1.5 mL	Nuclease-Free Water
Strip	24	8-channel strip
Protocol	1	Instruction guide for user
Base	2	A rack for 8 Auto Tubes

6. Auto Tube Content

Well	Buffer	Volume (µL)
1 / 7	Lysis Buffer	700
2 / 8	Washing Buffer 1	800
3 / 9	Magnetic Beads	800
4 / 10	Washing Buffer 2	800
5 / 11	Washing Buffer 2	800
6 / 12	Elution Buffer	100

7. Kit Storage and Shelf Life

- Components under room temperature (15~35°C) can be stored until the expiration date labeled on the box.
- The Proteinase K was transported at room temperature. When received, please store at 4°C.

8. Precautions

- For research use only.
- Avoid using expired reagents.
- When the temperature is below 20°C, place the Auto Plates / Auto Tubes in an oven (preheated 42~60°C) 5 to 10 minutes.
- Avoid vigorous shaking, in order to avoid excessive formation of foam.
- Carefully remove aluminum foil to avoid splashing.
- Do not expose the opened reagents or Auto Plates / Auto Tubes to air. The evaporation would lead to pH change, or effect on the extraction effectiveness.
- Please check the integrity of the Auto Plates / Auto Tubes and remember to insert the strips into the appropriate position of the suitable instrument before operating them.
- Please wear a mask and disposable gloves when handling.
- Use sterile consumables to avoid nuclease contamination.
- Reagent solution contains guanidine salt, avoid using bleach containing detergent.
- Avoid eyes, skin, and clothing contact with reagents. In case of any contact, flush with flowing water.

- If any serious incident occurs, please report to the manufacturer and the competent authority of the member state in which the user and / or the patient is established.

9. Materials required, Not Supplied

- TANBead® Nucleic Acid Extraction System
Model: SLA-16 / 32 / E13200 series (non-sterile)
- Disposable gloves
- Scissors, utility knives
- Micropipette, disposable tips (10 µL / 200 µL / 1000 µL)
- 1.5 mL microcentrifuge tube
- 15 mL / 50 mL conical tube

10. Sample Collection, Transportation, and Storage

■ Sample collection and storage

- Serum, whole blood
 - Serum specimens must be obtained from serum collection tubes, whole blood specimens must be obtained from sodium citrate or EDTA collection tubes.
- Specimen storage
 - Fresh whole blood specimens can be stored at room temperature for 6 hours.
 - After centrifugation, the serum sample can be stored at
 - Room temperature for 24 hours.
 - 2~8°C up to 7 days.
 - 20°C for long-term preservation.

■ Specimen transportation

Transportation of whole blood, serum specimens should be followed by specific pathogen transportation-related laws. The whole blood sample should be kept between 2~25°C during transportation and within 6 hours for separated serum. Serum samples can be transported between 2~8°C or by freezing.

11. Nucleic Acids Extraction Protocol

Before operating, turn on the warm-up system of TANBead® Nucleic Acid Extractor, if it is equipped with temperature controller, please setting at 45°C.

- Preparing the Assembled Auto Tubes by inserting Auto Tubes into the Base completely.
- Carefully remove the aluminum foil on the Auto Tubes.
- Pipet 200 µL serum or PBS suspension and 10 µL Proteinase K into column #1 / #7 of Auto Tube.
Note: The volume ratio of mixture and lysis buffer is about 200 µL : 700 µL. If it is changed, it might be affected the performance.
- Push Assembled Auto Tubes completely to the bottom of the plate rack. Make sure that the chamfer of the plate is at the lower left.
- Push strips completely to the bottom of strip rack frame.
- Close the door panel.
- Select the program "VIRUS-W4-AUTO". The parameters are given in following section.
- Carefully remove the Auto Tubes when the program is finished.
- Use micropipette to transfer the purified nucleic acid from column #6 / #12 to a clean tube.
- Discard used Auto Tubes and strips into the waste recycling bin.

12. Program

■ SLA-16 / 32 series

Program Name: VIRUS-W4-AUTO					Model: SLA-16 / 32 series			
Step	Well	Mixing (M)	Collect (S)	Rod	Mixing speed	Volume (μL)	Pause	Vapor (M)
1	3	1	60	On	Medium	800	Off	0
2	2	1	0	Off	Medium	800	Off	0
3	1	20	0	Off	Low	900	Off	0
4	2	0	60	On	Medium	800	Off	0
5	1	10	60	On	Medium	900	Off	0
6	2	2	60	On	Medium	800	Off	0
7	3	2	60	On	Medium	800	Off	0
8	4	2	60	On	Medium	800	Off	0
9	5	2	60	On	Medium	800	Off	10
10	6	5	120	On	Medium	150	Off	0
11	5	1	0	Off	Medium	800	Off	0
12	0	0	0	Off	Medium	0	Off	0

■ SLA-E13200 series

Program Name: VIRUS-W4-AUTO						Model: SLA-E13200 series			
Step	Well	Temp (°C)	Mixing (M)	Collect (S)	Rod	Mixing speed	Volume (mL)	Pause	Vapor (M)
1	3	45	1	60	On	Medium	800	Off	0
2	2	45	1	0	Off	Medium	800	Off	0
3	1	45	20	0	Off	Low	900	Off	0
4	2	45	0	60	On	Medium	800	Off	0
5	1	45	10	60	On	Medium	900	Off	0
6	2	45	2	60	On	Medium	800	Off	0
7	3	45	2	60	On	Medium	800	Off	0
8	4	45	2	60	On	Medium	800	Off	0
9	5	45	2	60	On	Medium	800	Off	10
10	6	45	5	120	On	Medium	150	Off	0
11	5	NA	1	0	Off	Medium	800	Off	0
12	0	NA	0	0	Off	Medium	0	Off	0

13. Result

Nucleic acid product purified by TANBead® nucleic acid extraction kit can perform qualitative / quantitative analysis of specific genes by PCR, RT-PCR, Q-PCR or Qrt-PCR. Please refer to the molecular diagnostic kit manual.

14. Reagent performance

■ Repeatability

A five-day reproducibility test was carried out with 100 IU / mL of EBV serum samples for 5 consecutive days with the same reagent kit by different operators. The coefficient of variation of nucleic acid extraction concentration is less than 5%.

■ Reproducibility

A five-day reproducibility test was carried out with 100 IU / mL of EBV serum samples for 5 consecutive days with the same reagent kit by different operators. The coefficient of variation of nucleic acid extraction concentration is less than 5%.

■ Detection limit of HCV virus: ≥100 IU / mL







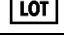
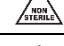





■ Interfering substance

The performance of this extraction kit will not be affected by EDTA, Li-Heparin, Sodium Citrate, D-Glucose, Hemoglobin, lipoprotein, and triglyceride in samples.

■ The stability of extracted RNA

Storage Conditions	RNA stability
-80°C	Over 90 days
-20°C	28 days
4°C	14 days
25°C	2 days
Freeze-thaw	5 times

15. Explanation of Symbols

	Manufacturer		Consult instructions for use
	Temperature limit		Contains sufficient for test
	Catalogue number		Caution
	Batch code		Non-sterile
	Do not re-use		Keep away from sunlight
	Date of manufacture		Use-by date
	For research use only		